



UNIVERSIDAD COMPLUTENSE DE MADRID
Facultad de Veterinaria
Departamento de Sanidad Animal



AFRICAN SWINE FEVER ANTIBODY DETECTION: IMMUNOBLOTTING ASSAY (IB)

jmvizcaino@vet.ucm.es
Av/ Puerta de Hierro s/n.
28040 Madrid.

Tel: (34) 913944082
Fax: (34) 913943908



1. MATERIALS

Watch video (Immunoblotting)

- Single channel pipettes 1-10 μ l
- Single channel pipettes 10-100 μ l
- Single channel pipettes 10-200 μ l
- Single channel pipettes 200-1000 μ l
- Chamber 37°C
- Analytical Balance
- Distilled water
- Minincubation trays (ref:170-3902.BIORAD)
- pH meter
- Tubes shaker or vortex mixer
- Reagent reservoir Polystyrene 50ml (ref: 4870.COSTAR)
- Sterile plastic tubes (10ml,50ml)
- Table centrifuge
- Latex or nitrile gloves

2. REAGENTS:

Watch video (Immunoblotting)

- **Blocking solution: 2% (w/v) milk in PBS/Tween-20 pH 7,2**

PBS Buffer pH 7.2/ 0.05%Tween-20:

ClNa [Merck 1.06404]-----	8 gr
ClK [Merck 1.04873]-----	0,2 gr
PO ₄ H ₂ K [Merck 1.06586]-----	0,2 gr
PO ₄ HNa ₂ [Merck 1.04936]-----	1,15gr
Tween-20 [Merck 8.22184]-----	0,5 ml
H ₂ O destilada -----	1000 ml

Check the pH before use. Store at 4°C

PBS/ Tween 20, 2% milk Buffer :

Non fat dry milk (NESTLÉ-Sveltesse o Molico)-----	2g
Buffer PBS pH 7.2 – Tween-20 0.05%-----	1000ml

Store at 4°C. Do not use it after two days.



➤ **Conjugate:**

Protein A peroxidase 1mg/ml (PIERCE, ref.0032400) store at 4°C. Resuspend it in 200 µl of distilled water and store it at -20°C. Before the addition to immunoblotting strips, dilute it 1/1000 in PBS 1x/0,05% Tween-20 /2% Milk.

➤ **Substrate solution:**

- a) Dissolve 6mg of 4-chloronaphtol in 2 ml of Metanol
- b) Add slowly 4-chloronaphtol/Metanol solution to 10 ml of PBS buffer pH 7.2 with vigorous agitation.
- c) Then, add 4 µl of H₂O₂ 30% (Panreac) to the PBS/4-chloronaphtol solution.

3. METHODOLOGY

Watch video (Immunoblotting)

3.1) Number the nitrocellulose strips.

3.2) Number the trays

3.3) Place the strips in the trays.

3.4) Block the nitrocellulose strips with blocking solution (1ml/strip) (PBS Buffer 7.2-Tween 20 0,05%-2% Milk)

3.5) Incubate for 30 minutes at 37°C in continuous agitation.

3.6) Wash four times with PBS buffer pH 7.2-Tween 20 0.05%-2% Milk, 1ml per strip.

3.7) Add 1ml of the test and control serum at a 1/40 dilution in blocking solution (PBS Buffer 7.2-Tween 20 0,05%-2% Milk).

3.8) Incubate it for 45 minutes at 37°C in continuous agitation.

3.9) Wash four times with PBS buffer pH 7.2-Tween 20 0.05%-2% Milk, 1ml per strip.



3.10) Preparation and addition of the conjugate:

*11µl Protein A + 10.989 11µl PBS Buffer pH 7.2-Tween-20 0,05%, 2% milk
(for un tray)*

3.11) Incubate it for 45 minutes at 37°C in continuous agitation.

3.12) Wash four times with PBS buffer pH 7.2-Tween 20 0.05%-2% Milk.

3.13) Substrate preparation:

6mg 4-chloronaphtol + 2ml Metanol + 10 ml PBS Buffer + 4 µl H₂O₂

3.14) 1ml of substrate is added

3.15) Stop the reaction after 10-15 minutes with distilled water.

4. READING AND INTERPRETING THE RESULTS

Sera showing a specific pattern of reaction similar as the antigen strips stained with positive control serum will be considered as positive to the ASF antibodies.

Any ELISA-positive or doubt serum that does not clearly react with the ASF proteins will be considered negative to ASF antibodies.

